Projectina AG Dammstrasse 2, P.O. Box CH-9435 Heerbrugg Schweiz/Switzerland Phone: +41-71-727 28 00 Fax: +41-71-727 28 28 marketing @projectina.ch www.projectina.ch



Cyanoacrylate fuming chamber with Touch Panel



The Cyanoacrylate fuming chamber, now available with touch panel, is a good method to secure evidence for all items with non-absorbent surface.

With the new touch panel and its simple operation, the evaporation process can now run as an automatic mode or continue to be operated manually.

In the automatic process, first the relative humidity will be generated up to 80%, and the process will begin signalled by an audible tone. During the actual vaporization process, it is important to manually watch for developments. After completion of the vaporization process, the air cleaning process inside the interior of the device can be activated by an easy click.

Health and safety during operation of the device are very important. The advanced security features will give you help, such as the automatic door lock from the beginning of the vaporization process, until the end of the air cleaning process. When to change the cyan filter blocks and fill the water tank will be displayed automatically.

Projectina AG

Dammstrasse 2, P.O. Box

CH-9435 Heerbrugg

FIUIIB. 347.7.22 28 28 marketing @projectina.ch Projectina AG Schweiz/Switzerland

Phone: +41-71-727 28 00 www.projectina.ch



Information about the fuming process:

The process of Cyanoacrylate fuming is a very efficient method to secure latent finger-prints on smooth, nonporous surfaces.

This method presents a considerable advantage compared to the securing of evidence with adhesions promoters, especially with older traces, and sticky or porous surfaces.

Cyanoacrylate (ethyl ester) is a viscid special adhesive that can be assigned to instant adhesives (super glue). Due to its ability of polymerisation, Cyanoacrylate is an extraordinary compact and fast hardening special adhesive. Fluid Cyanoacrylate generates vapours that react sensitively to certain constituents of latent finger-prints (sebum). This procedure is effective on nearly every surface.

With a two-chamber system and maintenance-free air purge cycle and no connection with an exhausting system necessary, simply connect to an electrical supply.

Technical data from the Cyanoacrylate fuming chamber

external dimension*:	
width in mm	900
height included the base in mm	2090
depth in mm	600

internal dimension*:	
width in mm	580
height in mm	1250
depth in mm	410

weight	approx. 159 kg
connected load	230 V, 1,3 Kw, 50 Hz
exhaust ventilator, type E 05	230 V / 140 W
- circulatory	10 m³/min
- total pressure difference	Pa 430
fan 2 pieces	230 V, 12 W
illumination	230 V, 36 W
vaporiser - heating plate	230 V, 350 W
humidification	230 V, 512 W
electronic temperature control unit	230 V
electronic moisture control unit	230 V
electronic humidity sensor	0-1 Volt
nominal current	6,0 A
fuse protection	16 A
control voltage	24 V DC
Intermediate shelf adjustable	
material	
- polypropylene cabinet	
- front door contains of laminated glass	
- hanging rods of stainless steel	

^{*}Various sizes available on request.

Made in Germany